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1. CONVENTIONAL USE

The equipment may only be used for the purpose as set forth in the catalogue and in the technical descriptions which means for ultrasonic cleaning as described in this instruction, and only in conjunction with external devices and components which are recommended or approved by EMAG AG. Any other use of the equipment is considered to be non conventional.

In case of non conventional use, EMAG AG will not be liable for possible damages and will not warrant for the correct and functional work of the components. Part of the conventional use is also the observance of all information given in the operating manual. To this attention is drawn by the  sign on the type plate. The correct and safe operation of the equipment implies appropriate transportation, storage, installation and assembly as well as careful handling and maintenance.

2. NON CONVENTIONAL USE

In case of non conventional use of the equipment danger for life and limb as well as danger for property may occur. Therefore a non conventional use is prohibited. In case of non conventional use EMAG AG will not be liable for possible damages and will not warrant for the correct and functionally work of the components. In order to avoid non conventional use of the ultrasound cleaning equipment the following directions are to be considered:

- The equipment may only be operated with the voltage printed on the type plate on the backside/ bottom side of the housing (shockproof socket).
- No flammable or aggressive chemicals (i.e. acids etc.) must be used for operation of the equipment. If necessary, use glass container EM 50 DG.
- During operation do not touch the liquid.
- In case of transportation damage, don't connect the equipment with main source.
- Do not clean creatures (i.e. finger nails, animals etc.) in the equipment.
- The equipment must only be operated with the correct level of water (water till mark, approx. 1 cm below upper brim). This must be controlled if equipment is operated for a longer time.
- Strongly contaminated goods shall be positioned in the equipment with the contaminated side to the bottom – do not staple goods.
- The goods to be cleaned must not be positioned directly onto the bottom of the tub – use the basket.
- If the equipment is operated for a longer time, do not stay longer than 4 hours within 1 m distance to the equipment.
- The equipment must only be operated indoor and within a dry environment.
- It is normal, that the upper brims warm up at longer operation.
- Before emptying the tub, the power connector must be disconnected from the main source.

3. OPERATING INSTRUCTION

Check the specification table for the correct operating requirements (located on the rear of the unit).

Startup

Before connecting power, put the unit on a stable and flat working surface. Add water to tank, suitable to completely immerse the items to be cleaned.

Connect the power cable on the rear of the unit and ensure the wall supply is switched „ON“. Touch „ON/OFF“ button on the front panel to turn on the machine. You are now ready to set the timer and temperature. Put the items to be cleaned in the basket. Switch on Ultrasonic with „ON/OFF“ button.

Setting the Temperature (optional at HC-Series, D-Series and H-Series)

When power is on, LED actual fluid temperature display window lamp flashes and displays fluid's actual temperature. To set related gears, rotate heating knob. In order to heat the cleaning fluid during cleaning, so long as the heating switch is rotated clockwise, then it will enter into the heating state.

The heating temperature can be selected and set freely according to the indication at the edge of rotary switch on the panel. Once the green heating working indicator lamp is on, the heating band PTC will also start working and heating the fluid. When the temperature reaches the set temperature, the green heating working indicator lamp will go off and the heating band PTC will stop working. When the fluid temperature is lower than the set temperature, the heating band PTC will also start working and heating the fluid. Then, the temperature of the fluid will be automatically controlled at the set temperature.

To stop heating, please rotate heating switch back to the „OFF“ position anticlockwise.

Important: Heating function is not restricted by the ultrasonic cleaning timing but can work for a long time (however, it is necessary to make sure there is enough water inside the slot).

Setting the timer

When power is on, rotate timing knob to set relevant gears. At this time, once „ON/OFF“ button is pressed, the ultrasonic cleaning function will start immediately. Then the LED window lamp flashes and displays the remaining time. When cleaning time is due, ultrasonic will stop and LED remaining time display window flashes and resumes to flash in the position before setting. At this time, as long as „ON/OFF“ button is pressed, cleaning will start again. This is the so-called “memory function”. It is convenient to use.

Ultrasonic function

After pressing „ON/OFF“ button each time, the machine will initiate Ultrasonic function and the ultrasonic will stop outputting. Press the button again to operate continuously after observing the cleaning effect.

Degassing function (optional at D-Series)

At the normal working state, once degas button is pressed, the machine will start degassing function and degas button indicator lamp will be on. Once degas button is pressed again, degassing function will stop and degassing Button indicator lamp will be off.

Sweep function (optional at D-Series)

At the normal work state, sweep function will start and sweep button indicator lamp will be on once sweep button is pressed. If sweep button is pressed again, the machine will stop sweep function and sweep button indicator lamp will be off.

Important: Sweep and degassing function do not work at once!

Emptying the tank

Remove waste liquids to empty the tank after no more need to use. Some models feature drain tap located on the side of the unit, using it can make the removal of the waste liquids effortless and safe. Always ensure the drain tap is “Closed” when filling the tank with water.

1. DESCRIPTION

**EMMI®-12HC, EMMI®-20H, EMMI®-20HC, EMMI®-30H, EMMI®-30HC, EMMI®-35HC-Q,
EMMI®-40H, EMMI®-40HC, EMMI®-55HC-Q, EMMI®-60H**

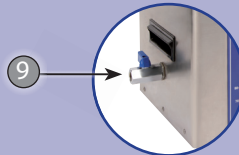
1. Cover – stainless steel
2. Basket - stainless
3. Tub – stainless steel
4. Power cable
5. ON/OFF
6. Timer
7. 3 stages ultrasound power
(Optional at Emmi®-12HC, Emmi®-20HC,
Emmi®-30HC, Emmi®-35HC-Q, Emmi®-40HC,
Emmi®-55HC-Q)
8. Switch for temperature setting regulator
9. outlet valve
(Optional at Emmi®-35HC-Q, Emmi®-40H,
Emmi®-40HC, Emmi®-55HC-Q, Emmi®-60H)

EMMI®-60HC, EMMI®-85HC, EMMI®-120HC, EMMI®-280HC, EMMI®-420HC

1. Cover – stainless steel
2. Basket - stainless
3. Tub – stainless steel
4. Power cable
5. ON/OFF
6. Timer
7. ultrasound power regulator
8. Switch for temperature setting
9. outlet valve

2. TECHNICAL DATA

All units have a power supply of 230V ~ / 50Hz and the protection class IP20.

The heating controller has variable control in the range between 20 - 80 ° C.

	External dimensions LxWxH (mm)	Internal dimensions LxWxH (mm)	Content (liter)	Ultrasound power (watt)	heating power (watt)	Generator	outlet valve
Emmi [®] -12HC	225 x 125 x 170	200 x 100 x 70	1,2	80	200	2	no
Emmi [®] -20H	260 x 150 x 210	230 x 118 x 80	2,0	150	200	2	no
Emmi [®] -20HC	260 x 150 x 220	230 x 120 x 80	2,0	150	200	2	no
Emmi [®] -30H	265 x 160 x 230	240 x 135 x 100	3,0	200	300	3	no
Emmi [®] -30HC	265 x 165 x 240	240 x 135 x 100	3,0	200	300	3	no
Emmi [®] -35HC-Q	325 x 260 x 210	300 x 240 x 65	3,5	250	150	6	yes
Emmi [®] -40H	325 x 175 x 230	300 x 155 x 100	4,0	250	300	4	yes
Emmi [®] -40HC	325 x 180 x 140	300 x 155 x 100	4,0	250	300	4	yes
Emmi [®] -55HC-Q	325 x 260 x 290	300 x 240 x 150	5,5	300	300	6	yes
Emmi [®] -60H	235 x 175 x 285	300 x 155 x 130	6,0	350	450	6	yes
Emmi [®] -60HC	325 x 175 x 285	300 x 155 x 130	6,0	350	450	6	yes
Emmi [®] -85HC	315 x 250 x 250	300 x 240 x 150	8,5	400	400	4	yes
Emmi [®] -120HC	315 x 250 x 300	300 x 240 x 200	12,0	800	450	6	yes
Emmi [®] -280HC	530 x 320 x 360	500 x 300 x 200	28,0	1000	1000	8	yes
Emmi [®] -420HC	540 x 340 x 400	500 x 300 x 300	42,0	1500	1350	12	yes

Technical changes reserved.

4. DESCRIPTION

D-SERIE

EMMI®-D20Q, EMMI®-D30, EMMI®-D40, EMMI®-D60, EMMI®-D100, EMMI®-D130, EMMI®-D280

1. basket - stainless
2. tub – stainless steel
3. power cable
4. heating
5. timer
6. 50% Power
7. sweep
8. degas
9. start/stop
10. on/off

without image: Cover – stainless steel
11. outlet valve

H-SERIE

EMMI®-05ST, EMMI®-08ST, EMMI®-08STH, EMMI®-H12Q, EMMI®-H20Q, EMMI®-H22, EMMI®-H30, EMMI®-H40, EMMI®-H60, EMMI®-120L, EMMI®-H420, EMMI®-H280

1. basket - stainless
2. tub – stainless steel
3. on/off
4. heating-control lamp
5. heating
6. timer
7. ultrasound-control lamp

without image: Cover – stainless steel
8. outlet valve

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6. EXAMPLES FOR THE CLEANING OF GOODS WITH ULTRASOUND

- Eyeglasses, glass frames (please pay attention: Glasses must not be in direct contact with the tub bottom).
- Gold and silver jewels, jewelry, emeralds, pearls
- Ink crayons, drawing instruments / -accessories
- Dentures
- Aircraft components, engine components, electrical boards, ink cartridges, motor bike chains, tools, carburators, etc.
- Precision parts, weapon parts, i.e. magazines, bullet casings, strikers, springs.

7. CONCENTRATE FOR DISOXIDATION

For the cleaning operations various cleaning concentrates are available which can make cleaning more efficient. These are available at the place where you have bought your equipment or through www.emag-germany.de

EMAG recommends for the various requirements the following high efficient concentrates:

- EM 070** for dental application and cleaning of dentures at home. Application: 1%.
- EM 080** Eyeglasses, jewelry, optical glasses, ink crayons smaller tools and instruments such as small precision parts, shaver heads. Application: 1%. (except for contact lenses).
- EM 100** Concentrate for disoxidation of silver etc.). Application: neat.
- EM 200** Concentrate for disinfection of instruments. Application: 1%.
- EM 300** Super plus – extreme cleaner for difficult cases (supports, hinges, weapon parts, bullets, brass etc.). Application: 1%.
- EM 303** PC Board cleaning. Application: 5-10%.
- EM 404** Aluminum and diecast cleaning (carburetors, engine blocks, cylinder heads). Application: 3-5%.
- EM 600** Spezialreiniger (Glas, Gummi, Keramik, Kunststoff, Metall). Application: 1-10%.
- EM 700** Non-ferrous metal cleaner (Bullet casings, PC boards from maintenance, Jewellery, Clock parts). Application: 5%.

If concentrates are used, application instructions must be followed very correctly!

If cleaning concentrates of other manufacturers are used, we will not be liable for any problems occurring with the goods to be cleaned and any damages which may happen to the equipment. For detailed information you can contact us at any time.

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8. WARRANTY

This product is warranted against defects in manufacturing or materials for a period of two year from the date of purchase. Defects may be corrected by repair, or by replacement of the unit, at the option of the manufacturer.

This warranty may be voided if the unit is not operated in accordance with the instructions provided by the manufacturer. The warranty may also be voided by the opening of the unit, and/or its partial or total assembly, without prior factory authorization.

Any claim made under this clause shall be directed to the seller, and shall be verified by proof of purchase. The seller will advise the buyer concerning disposition of the unit. The normal procedure shall be for the manufacturer to issue a warranty claim number, and to authorize return of the unit to the factory.

Buyer shall bear shipping costs.

© Copyright EMAG AG. All rights reserved. Manufactured in the European Community. The content of this Operating Instruction is subject to the copyright and it must not be copied neither in part nor in total. Changes, reductions, extensions and amendments of any publication or translation requires the prior written agreement of EMAG AG.

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© Copyright EMAG AG. Alle Rechte vorbehalten. Hergestellt in der BRD/EU. Der Inhalt dieser Bedienungsanleitung unterliegt dem Urheberrecht und sie darf weder ganz noch teilweise kopiert werden. Veränderungen, Kürzungen, Erweiterungen und Ergänzungen jede Veröffentlichung oder Übersetzung durch Dritte bedarf der vorherigen schriftlichen Einwilligung durch EMAG AG.